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Warranty

In accordance with IKA® warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs. The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

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The tool kit is part of the delivery scope of the disperser.

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Assembling the dispersion tool (overview Fig. 2)

- Open the knurled screw (item 3) so that the thread does not protrude into the hole.
- Insert the dispersing element (item 2) into the drive unit (item 1) as far as the stop. After some resistance (depending on the model of the drive - pressure ball) you will hear the shaft clicking into place. The dispersing element is fitted correctly if the recessed edge of the dispersing element is flush with the bottom edge of the drive unit.

- **Attention:** Now secure the dispersing element by screwing the knurled screw tight (item 3).

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IKA Japan K.K.
Phone: +81 6 6730 6781
eMail: info_japan@ika.ne.jp

INDIA

IKA India Private Limited
Phone: +91 80 26253 900
eMail: info@ika.in

UNITED KINGDOM

IKA England LTD.
Phone: +44 1865 986 162
eMail: sales.england@ika.com

VIETNAM

IKA Vietnam Company Limited
Phone: +84 28 38202142
eMail: sales.lab-vietnam@ika.com

Thailand

IKA Works (Thailand) Co. Ltd.
Phone: +66 86 375 7451
eMail: sales.lab-thailand@ika.com

TURKEY

IKA Turkey A.Ş.
Phone: +90 216 394 43 43
eMail: sales.turkey@ika.com

Discover and order the fascinating products of IKA online:
www.ika.com

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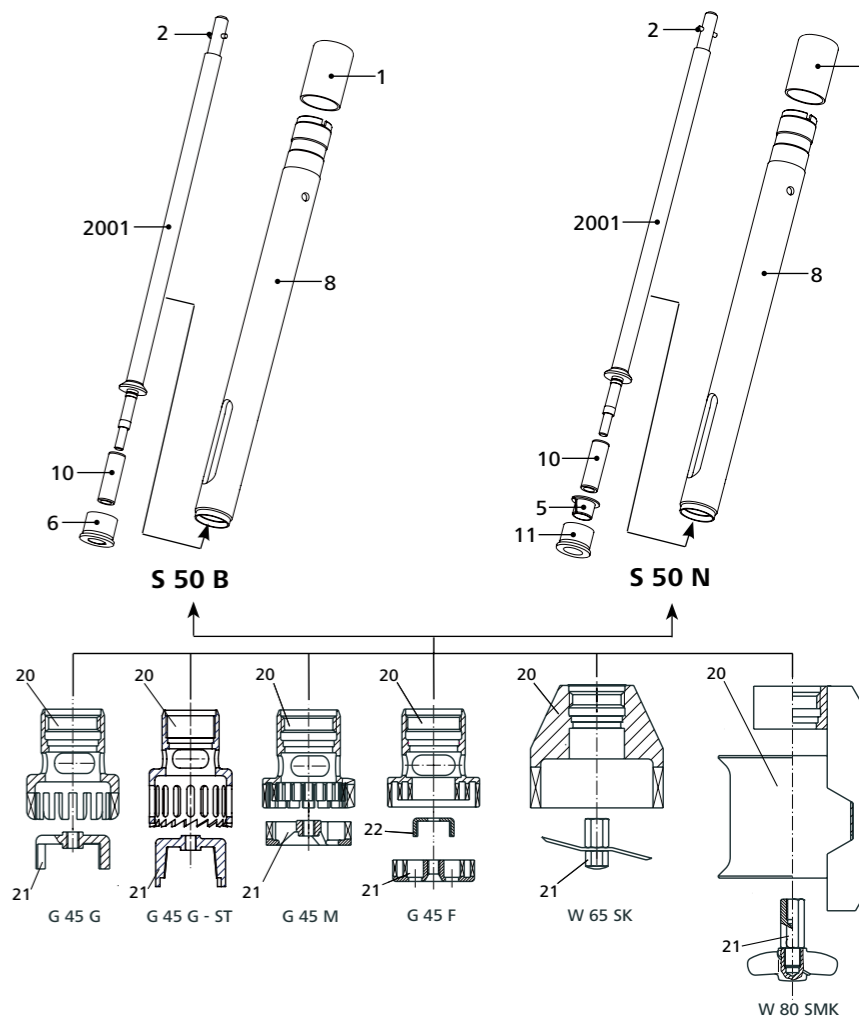
Safety instructions EN

- Read the operating instructions in full before starting up and follow the safety instructions.**
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

- ⚠

The dispersion tool may heats up during operation. In particular in the processing of hot medium, the tool must be cooled before it can be removed from the drive.
- When in operation the dispersion tool can become hot. If the dispersion tool is not inserted into the drive flange correctly or the medium is hot, it can become extremely hot.

- Wear your personal protective equipment in accordance with the hazard category of the medium to be processed. Otherwise there is a risk of splashing and evaporation of liquids; body parts, hair, clothing and jewelry getting caught.
- The work has to be interrupted immediately if you notice unusual noise and/ or increasing emission of liquid from the upper, sidewise rinsing drilling. The bearings must be controlled, and replaced if necessary.
- Note the operating instructions of the disperser unit.
- Abrasion of the dispersion equipment or the rotating accessories can get into the medium you are working on.
- Do not use any damaged dispersion tools e.g. with hairline cracks or splits.
- Attention!** Never run the dispersion tool dry. Without cooling by the medium the bearing will be destroyed.
- The dispersion tool is not suitable for continuous operation. Likewise the optimal dispersion duration and rotating frequency must be determined by attempts. Usually a few seconds are sufficient, maximum duration is one minute. Longer application times bring no improvements, increase however the sample temperature substantially.
- Dispersing causes the medium to heat up.
- The fill level of the medium must be at least 20 mm below the upper hole of the shaft.
- The function of the dispersion tools depends on the condition of the sharp edges of rotor and stator (integrated in the shaft tube). Abrasive media can round off these edges fast, whereby the disperse effect diminishes.
- For **S 50 N**, the gasket and bearings are made of PTFE and stainless steel; the following points should therefore be noted: *Chemical reactions of PTFE occur in contact with molten or solute alkali metals and alkaline earth metals, as well as with fine powders of metals in groups 2 and 3 of the periodic system at temperatures above 300 °C - 400 °C. Only elementary fluorine, chlorotrifluoride and alkali metals attack it; halogenated hydrocarbons have a reversible swelling effect.* (Source: Römpps Chemie-Lexikon and "Ulmann", Volume 19)

Product information

- The shaft bearing of the dispersion tool consists of a slide bearing (**S 50 N** – PTFE, **S 50 B** – bronze).
- Content of package: Shaft and dispersion head.
- The generator (dispersion head) consists of a rotor and a stator. These parts can be unscrewed from the shaft. Note that rotors and stators form matching pairs and must not be confused. Five different dispersion heads are available for various applications.
- The materials used conform to the FDA (Food and Drug Administration Approval) - rules.

Application instructions

Application: Wet crushing, suspension, emulsifying (Batch-operation).

Spare parts list

Item	Designation	Item	Designation	Item	Designation	Spare parts diagram see front page.
1	Plastic cap	8	Shaft	21	Rotor	For orders of spare parts, please specify the type of dispersion tool and the designation of the spare part.
2	Grooved pin	10	Shaft sleeve	22	Blade	
5	Bearing bushing (PTFE)	11	Shaft sleeve	2001	Rotary shaft	
6	Bearing bushing (bronze)	20	Stator			

Cleaning, disinfection and sterilization

Allowed procedures	Sterilization techniques
Damp heat	Autoclave to 121 °C at 2 bar positive pressure.
Chemical procedures	By germ - killing dissolvent alcohol, phenol, formalin..., disinfectant remainders are to be removed with germ - free water.
Hot air	Germ - killing by hot-air with 160 to 190 °C (approx. 30 min.).

For pre-cleaning the dispersion tool, operate it in a detergent for dissolving the residuary of the substance. Due to the large flow rate, rotor and stator are cleaned to a large extent. The dispersion tool must be divided and cleaned immediately after working, so that the adhering residuary of the substance don't cultivate unwanted bacterial cultures.

For cleaning purposes disassemble the dispersion tool as described below by the shaft and the generator key, both included in the tool kit of the dispersion drive:

- Unscrew the rotor (21) and blade (22) from the rotary shaft (2001) using the generator and shaft wrenches included in the tool kit. The generator wrench is used to hold the rotor while unscrewing it by turning the shaft from above with the shaft wrench.
- Caution:** The dispersion head **G 45 G** has two prominent blades.
- Unscrew the stator (20) from the shaft (8) (left-hand thread). Insert the cross pin (3 mm) of the shaft wrench into the slot on the end face of the shaft tube. The stator can now be unscrewed by turning it in the clockwise direction with the generator wrench.
- Pull the rotary shaft (2001) down and out of the shaft tube. The components (5, 6, 10 and 11) threaded onto the shaft can simply be pulled off.
- The bearing bushing (5 and 6) must be replaced if it is worn.

Reassemble the dispersion tool in the reverse sequence. When assembling ensure that the bearing bushing and the rotor are correctly installed. Attention! Do not tighten the rotor too strongly, because otherwise the small teeth at the rotor become bent and the disper­sion tool will be destroyed.

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