IKA

designed for scientists

R 472 Floor stand
R 474 Telescopic stand
T 653 Telescopic stand

Betriebsanleitung	DE	3
Operating instructions	EN	6
Mode d'emploi	FR	9

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Contents

	Page
EU Declaration of conformity	2
Explication of warning symbols	7
Safety instructions	7
Unpacking	8
Correct use	8
Set-up and Commissioning	8
Maintenance and cleaning	9
Accessories	9
Technical data	10
Warranty	10

Explication of warning symbols

Indicates an (extreme) hazardous situation, which, if not avoided, will result in death, serious injury.

Indicates a hazardous situation, which, if not avoided, can result in death, serious injury.

Indicates a potentially hazardous situation, which, if not avoided, can result in injury.

Indicates practices which, if not avoided, can result in equipment damage.

Safety instructions

- **Read the operating instructions in full before starting up and follow the safety instructions.**
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the device.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Set up the stand in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.

- **R 474 / T 653:** The stand pillar must be inserted into the baseplate so that the guide groove is at the back. Other alignments are not permissible, as otherwise the stand can topple over when a unit is mounted on it.

- **R 474 / T 653:** Any small unevennesses that cause the stand to rock must be compensated for with the height-adjustable feet.
- **R 472:** The castors must be locked when the device is in use.

- The **T 65 basic / digital ULTRA-TURRAX®** dispersion unit must be lifted on to the guide head of the telescopic stand with the aid of an assistant, and secured with the two screws and toothed lock washers (supplied), in accordance with the **T 65 basic / digital ULTRA-TURRAX®** operating instructions. The unit must be mounted in front, vertically above the baseplate; a different alignment is not permissible, as otherwise the assembly could topple over.
- The overhead stirrer **RW 47 digital** must be lifted on to the guide head of the telescopic stand with the aid of an assistant, and secured with the four screws (supplied), in accordance with the **RW 47 digital** operating instructions. The unit must be mounted in front, vertically above the baseplate; a different alignment is not permissible, as otherwise the assembly could topple over.

- This telescopic stand must be fitted with a safety disconnection device, which stops the unit if the stand is moved upwards when the unit is operating.
- Switch the unit off before moving the telescopic stand upwards.
- When lowering the telescopic stand, take care that no parts of the body get between the safety switches (danger of crushing).
- Only switch the unit on when the tool (agitator or dispersion unit) is in the vessel.

- The handwheel bolt on the stand can come loose due to vibration, which would lead to the unit moving down. So, check the tightness of the handwheel bolt regularly when using the unit, and retighten it if necessary.
- Raise or lower the stand (after slackening the handwheel bolt) by applying the manual force as close as possible to the stand pillar. The further the force is applied from the pillar, the more force you need to apply, plus there is an increased risk of the equipment toppling over.

NOTICE

- Secure the vessel against spinning round, by using a clamp, for example.
- Observe the operating instructions of the **T 65 basic / digital ULTRA-TURRAX®** dispersing unit or the **RW 47 digital** stirring unit!
- Check the appliance and accessories beforehand for damage each time you use them. Do not use damaged components.
- Protect the appliance and accessories from bumps and impacts.

Unpacking

Unpacking:

- Please unpack the device carefully.
- In the case of any damage a detailed report must be sent immediately (post, rail or forwarder).

Delivery scope:

Telescopic stand

- **R 472 Floor stand** or **R 474 Telescopic stand** or **T 653 Telescopic stand** according to order
- Operating instructions.

For technical reasons associated with shipping, the telescopic stand is supplied stripped down, as pre-assembled modules.

Intended use

Use:

The **R 472 Floor stand / R 474 Telescopic stand** is used solely for mounting **RW 47 digital** agitator units.

The safety of you cannot be guaranteed:

- If the device is operated with accessories that are not supplied or recommended by **IKA**.
- If the device is operated improperly or in contrary to the **IKA** specifications.

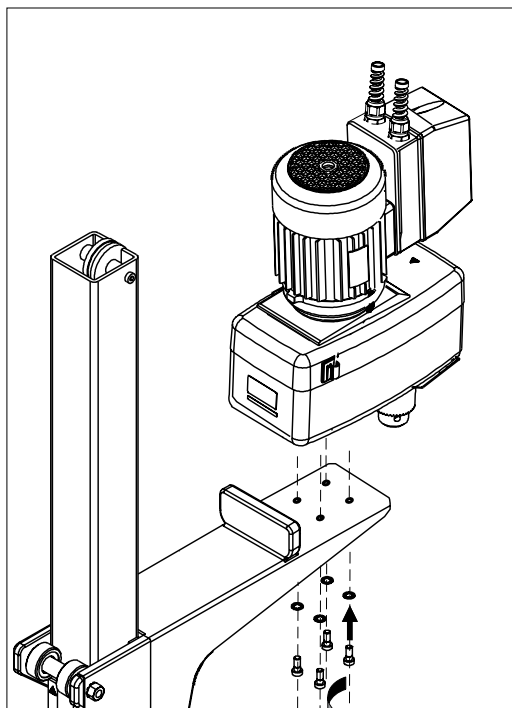
WARNING

R 472 Floor stand: the castors must be locked when the device is in use.

The **T 653 Telescopic stand** is used solely for mounting **T 65 basic / digital ULTRA-TURRAX®** dispersion units.

Set-up and commissioning


Setting up the stand (R 472 Floor stand):

Place **RW 47 digital** overhead stirrer on the arm of the stand. Fasten the overhead stirrer to the stand from below with hexagonal socket screws supplied. The help of an assistant is needed for screwing on the overhead stirrer. The cable clips included in the delivery package are required for fixing the power cable of the unit.

Check that the stirrer is held in position firmly prior to each use and also at regular intervals. The position of the stirrer must only be adjusted when the equipment is stationary and the power supply is disconnected.

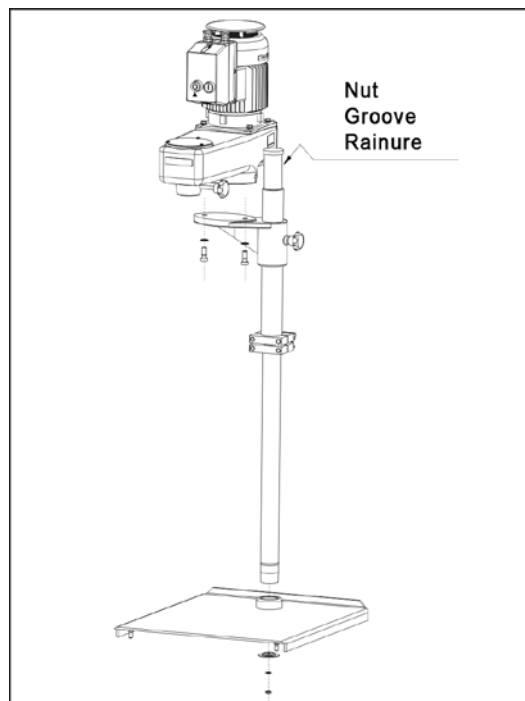Setting up the stand (R 474 / T 653 Telescopic stand):

The stand pillar is inserted into the baseplate so that the guide groove is at the back. The black handwheel bolt is then on the right. The stand pillar is then secured from below with a hexagon nut, toothed lock washer and a large shaped washer.

All four rubber feet are height-adjustable, so the stand can be set up so it does not rock, even on uneven floors.

In addition, a hole is provided in the baseplate of the **R 474 Telescopic stand**, for screwing on a 16 mm diameter support rod. A clamp can be fixed on this rod.

Mounting the unit:

The **RW 47 digital** agitator unit or the **T 65 basic / digital ULTRA-TURRAX®** dispersion unit are placed on the arm of the stand and bolted on from below with the bolts supplied. The help of an assistant is needed for screwing on the **T 65 basic / digital ULTRA-TURRAX®** dispersion unit.

The cable clips included in the delivery package are required for fixing the power cable of the unit.

Safety disconnection device:

The height-adjustable telescopic stand must be fitted with a safety disconnection device. We recommend installation of the **SI 400** safety limit switch together with the **SI 472 / SI 474** clamp support. You will find information on their installation in the operating instructions of the units.

Commissioning (R 474 / T 653 Telescopic stand):

The stand can be raised and lowered by loosening the handwheel bolt. The integral gas shock absorber helps to keep the manual force needed for this as light as possible. The manual force must be applied close to the stand pillar, so that the stand can be raised and lowered with only a slight use of force. If the force for raising or lowering is applied at a distance from the stand pillar, a turning moment results, leading to increased friction on the pillar guide. As a result, the stand can only be moved with great difficulty, or cannot be moved at all. Besides this, the danger of the stand toppling over is increased.

When the unit is at the correct height, the handwheel bolt is tightened by hand. Check the tightness of the handwheel bolt regularly when using the unit, and retighten it if necessary.

Maintenance and Cleaning

R 474 / T 653 Telescopic stand:

The stand pillar must always be lightly oiled, so that raising and lowering is not made more difficult due to increased friction.

Apart from this, the telescopic stand is maintenance-free.

Cleaning:

- Use only cleaning agents which have been approved by IKA to clean IKA devices: these are water (with tenside) and isoproApanol.
- Wear protective gloves during cleaning the devices.
- Before using another than the recommended method for cleaning or decontamination, the user must ascertain with IKA that this method does not destroy the device.

Spare parts order:

When ordering spare parts, please give:

- device type.
- serial number, see type plate.
- item and designation of the spare part, see www.ika.com.

Repair:

Please send the device for repair only after it has been cleaned and is free from any materials which may constitute a health hazard.

For repair, please request the "**Decontamination Certificate**" from IKA, or download printout of it from the IKA website www.ika.com. If you require servicing, return the device in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

Accessories

SI 400	Safety limit switch
SI 472	Support holder
SI 474	Support holder

Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

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IKA India Private Limited
IKA England LTD.
IKA Vietnam Company Limited