

# Thermometer Manual

## ST-10

- I. General Description  
Pen Thermometer which has the advantages of being multi-function, low power, high stability and high accuracy. It is a simple and practical high-temperature testing instrument.
  
- II. Product Functions  
Temperature measurement  
Range:  $-50^{\circ}\text{C} \sim +300^{\circ}\text{C}$  ( $058^{\circ}\text{F} \sim +572^{\circ}\text{F}$ )  
Temperature Accuracy:  $\pm 1^{\circ}\text{C} / 1.8^{\circ}\text{F}$   
Power Supply: 1.5V (AG13)\*1.
  
- III. Description of Button Functions
  - a. ON/OFF Button: Switch Function
  - b.  $^{\circ}\text{C} / ^{\circ}\text{F}$  Button: When the LCD displays a temperature, press to change display from  $^{\circ}\text{C}$  to  $^{\circ}\text{F}$
  - c. HOLD Button: When the LCD displays a temperature, press this button to hold the measured temperature on the display. Pressing the button again will revert the device back to temperature measurement. When implemented the hold symbol will be displayed on the LCD screen.