

EnSURE™ Touch

Simplifying Food Safety




Introducing EnSURE™ Touch

EnSURE™ Touch is the next generation ATP and environmental monitoring system. Designed to work like your smartphone, it effortlessly provides you with rapid and accurate data to support your food safety program. EnSURE Touch adapts to your workplace and is built with you in mind. Whether you're new to food safety or if you've been here a while, EnSURE Touch helps make everything a little bit easier.

How it works.


EnSURE™ Touch is a smart handheld instrument that rapidly measures contamination from bacteria, pathogens, foods, enzymes and indicator organisms. After taking a sample, the system measures the amount of Adenosine Triphosphate (ATP – found in all living or once living things) and gives you an objective numerical value combined with a Pass/Caution/Fail notification.

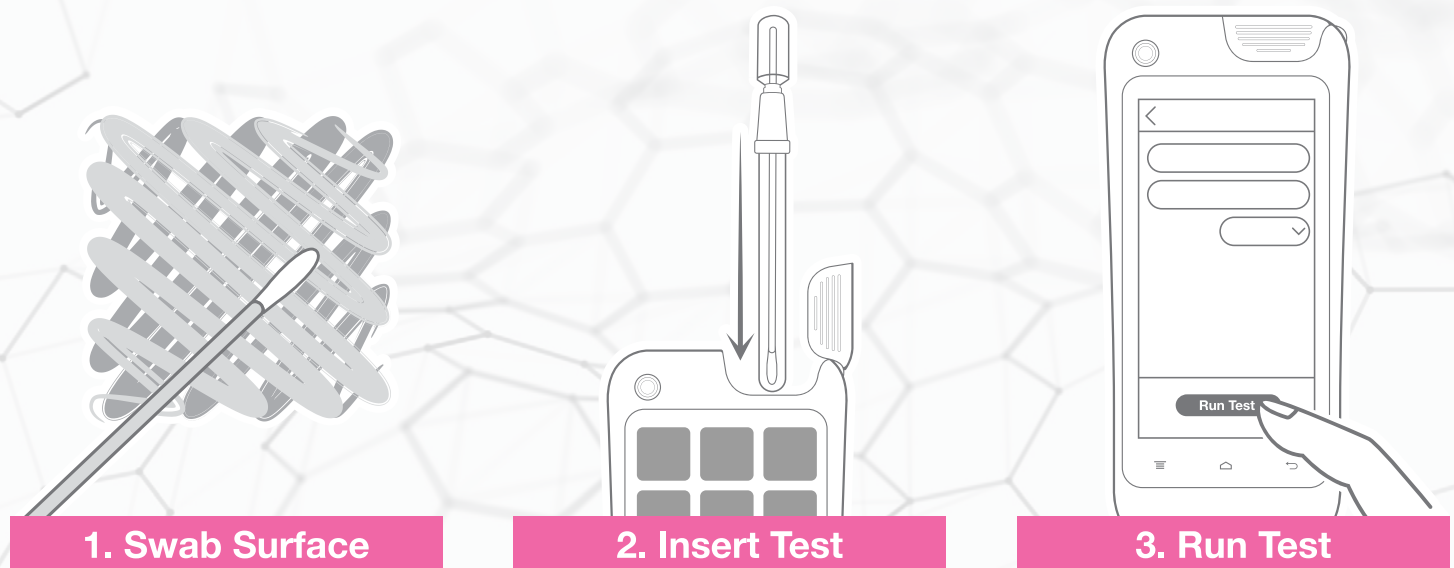

How can it help?

- Verify sanitation in 10 seconds
- Keep food safety data in one place
- Make audits easier
- Compare the effectiveness of sanitation procedures
- Objectively prove standards have been met
- Communicate success with insightful reports
- Identify problem areas
- Provide on-the-spot training and feedback
- Meet FSMA, GFSI, and ISO standards
- Protect brand reputation
- Reduce risk of food poisoning, facility closure, and product recalls
- Lower food safety costs by bringing testing in-house
- Provide evidence of due diligence – HACCP
- Optimize processes and limit down-time
- Makes environmental monitoring as easy as using a smartphone

Innovation at Your Fingertips

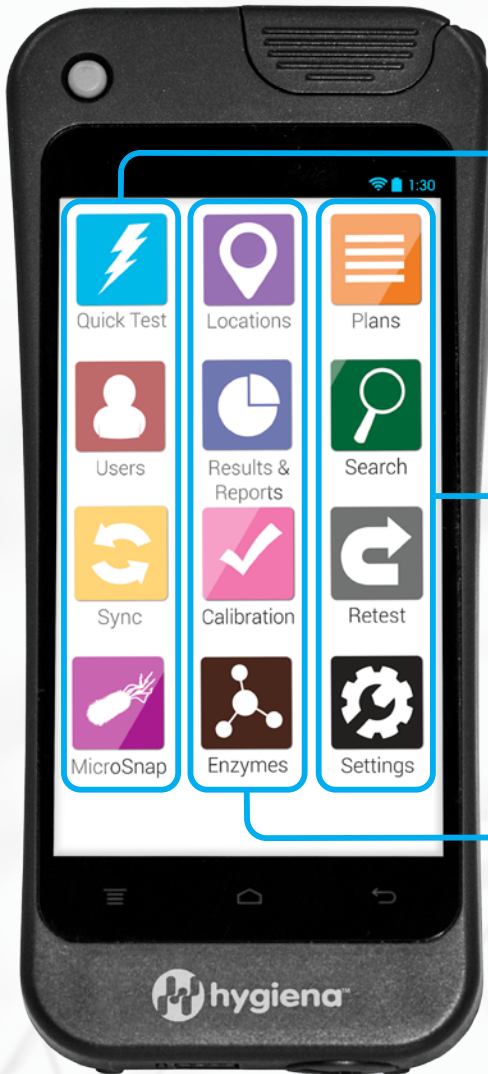

Quick Test
Run a test in two touches.



Sync
Wirelessly sync data to SureTrend Cloud.



Plans
Easily group, schedule, and randomize testing locations to ensure sampling coverage.



Retest
Swipe through failed tests and instantly run re-tests to show effectiveness of corrective action.



Locations
Create & edit test points with custom fields and pass/fail limits on the go.



Calibration
Verify calibration in just a few taps without having to send back to manufacturer.



Users
Manage access with password protection and roles.



MicroSnap
Test and show results for Coliforms, *E. coli*, *Enterobacteriaceae*, and Total Viable Count.



Search
Quickly find the location or plan you're looking for.



Settings
Customize settings for language, custom fields, and test types.



Results & Reports
Easily view testing results and trend analysis.



Enzymes
Test and show results for Alkaline Phosphatase and Acid Phosphatase.



The AOAC Research Institute, one of the world's most recognizable product certification organizations, issued its *Performance Tested Method*™ certificate to UltraSnap™ Surface ATP Tests for use with the EnSURE™ Touch.

Specifications:

- Size:** 7" x 3" x 1"
- Weight:** 0.7 lbs
- Sensitivity:** <1 femtomole of ATP
- Memory:** 2GB internal + cloud storage
- Connectivity:** USB or Wi-Fi
- Language:** Supports multiple languages
- Battery:** Rechargeable Lithium-Ion battery with USB-C charging*

* Battery life varies based on use and configuration

Always Up-To-Date Always Backed Up Always Encrypted




SureTrend Cloud is an easy-to-use online data analysis platform that syncs wirelessly with EnSURE Touch to unite testing data and bring it to life. It enables users to monitor testing results and manage user access across multiple facilities, making risk management easier than ever.



Features & Benefits:

- Easy-to-use & intuitive – no training required
- Web-based platform – no software downloads needed
- Unite multiple types of food safety data in one place
- Invite colleagues and assign roles to increase collaboration, accountability, and organizational buy-in and provide greater visibility into food safety program
- Interactive dashboard gives a birds-eye-view of overall testing performance
- Ready to use reports save time, custom reports add extra detail & personalization
- New capabilities added regularly via seamless and automatic software updates – no need to call IT
- Unlimited data and storage – complimentary startup costs
- Easily manage one or hundreds of instruments from one account
- Track re-tests to discover opportunities for retraining or identify equipment that may need replacing

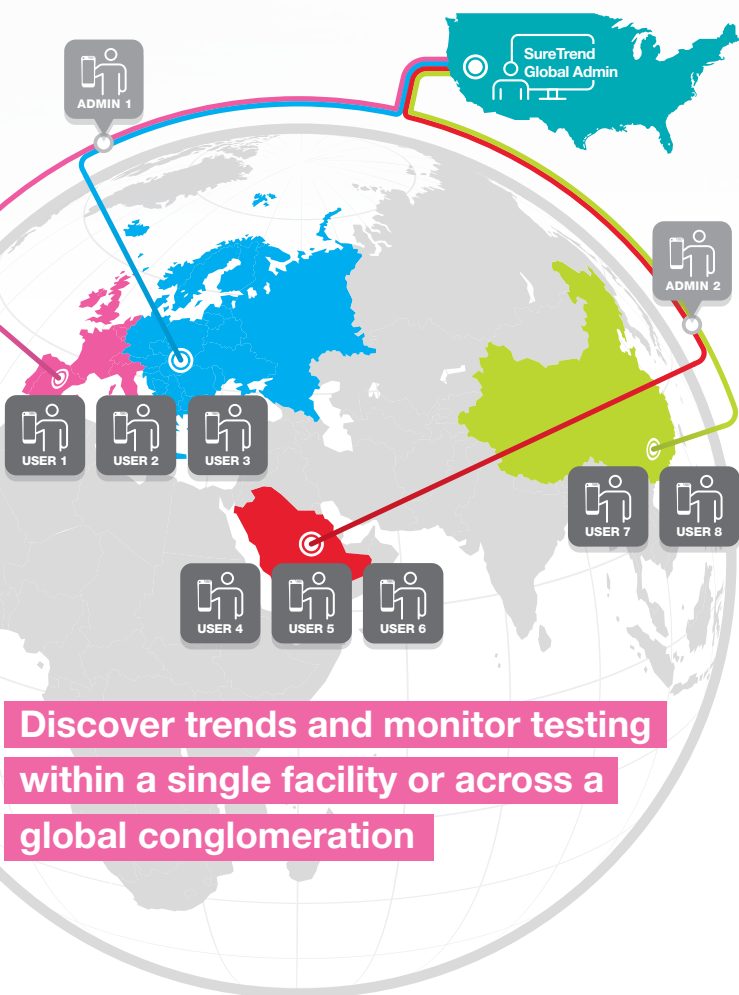

Discover trends and monitor testing within a single facility or across a global conglomeration

Test Devices

Hygiene test devices are unbelievably easy to use. Their design maximizes sample collection and recovery while using patented Snap-Valve™ technology to make activation easy. These devices eliminate extraneous materials, improve efficiency, minimize variability, and lower costs.

UltraSnap™ | Surface ATP Test



Catalog No. US2020 | Qty: 100

SuperSnap™ | High-Sensitivity Surface ATP Test



Catalog No. SUS3000 | Qty: 100
Catalog No. SUS3000X | Qty: 100

AquaSnap™ TOTAL | Water ATP Test



Catalog No. AQ-100X | Qty: 100

AquaSnap™ FREE | Water ATP Test



Catalog No. AQ-100FX | Qty: 100

CalCheck | LED Calibration Verification Device



Catalog No. CAL | Qty: 1

MicroSnap™ | Indicator Organism Tests



Catalog No. MS1-CEC | Qty: 100

Catalog No. MS2-ECOLI | Qty: 100

Catalog No. MS2-COLIFORM | Qty: 100

Catalog No. MS1-TOTAL | Qty: 100

Catalog No. MS2-TOTAL | Qty: 100

Catalog No. MS1-EB | Qty: 100

Catalog No. MS2-EB | Qty: 100

ZymoSnap™ ALP | ALP Pasteurization Verification Test



Catalog No. ZS-ALP-100 | Qty: 100

CROSS CHECK | ACP Cooking Efficiency Test



Catalog No. CX-3000 | Qty: 100

EnSURE Touch is factory calibrated and uses advanced technology that does not require recalibration.

CalCheck is used for in-house calibration checks as part of a quality control program so you don't need to mail it in for an annual calibration check.



MPB-049 REV D

Americas

Hygiena Headquarters

941 Avenida Acaso Camarillo, CA 93012
Phone: 1-805-388-8007
Fax: 1-805-388-5531
info@hygiena.com

Hygiena Canada

2650 Meadowvale Blvd Unit 14
Mississauga, Ontario L5N 6M5
Phone: 1-833-494-4362 (Toll-free)
or 1-416-686-7962

Hygiena Mexico, S.A. de C.V.

Calle 3 Anegas 409 Bodega 5
Col. Nueva Industrial Vallejo
Delegación Gustavo A. Madero
C.P. 07700, CDMX, México
Phone: +52-5281-4108 or +52-5281-4146

International

Hygiena International

8, Woodshots Meadow
Watford, Hertfordshire
WD18 8YU, UK
Phone: +44 (0)1923-818821
Fax: +44 (0)1923-818825

Hygiena (Shanghai) Trading Co., Ltd.

Rm.7K, No.518, Shangcheng Rd.
Pudong New District
Shanghai, China 200120
Phone: +86 21-5132-1081, +86 21-5132-1077,
+86 21-5132-1078
Fax: +86 21-5132-1080

Hygiena Diagnóstica España S.L.

P. I. Parque Plata,
Calle Cañada Real 31-35
41900, Camas, Sevilla, Spain
Phone: +34 954-08-1276

help.hygiena.com

 youtube.com/hygienatv

 vimeo.com/hygienatv

Hygiena™, EnSURE™, UltraSnap™, SuperSnap™, AquaSnap™, and MicroSnap™, are trademarks of Hygiena LLC.

